

TestSure 



PREMIUM UNIVERSAL

PH TEST STRIPS 0-14

EASY DIRECTIONS FOR USE

Grip A Single Test Strip From The Container, Hold The End Away From Colored Pads. Carefully Dip The Strip Into Any Liquid Or Gel For Half-Second. Allow All Four Pads To Get Wet And Then Remove From The Solution - Colors Will Change Almost Immediately. Match Panels To The Ph Diagram On The Box To Get Ph Reading. Ph Ranges From 0 (Very Acidic) To 14 (Very Alkaline). Neutral Is 7.

IMPORTANT:

- ◆ Keep away from light and moisture.
- ◆ Promptly replace cap after taking out strips.
- ◆ Do not remove desiccants.
- ◆ Do not touch test areas of reagent strips.
- ◆ Use it within the expiration date.
- ◆ Read insert carefully before use.

